


# Clean Test

Rapid test for the hygiene control of work surfaces

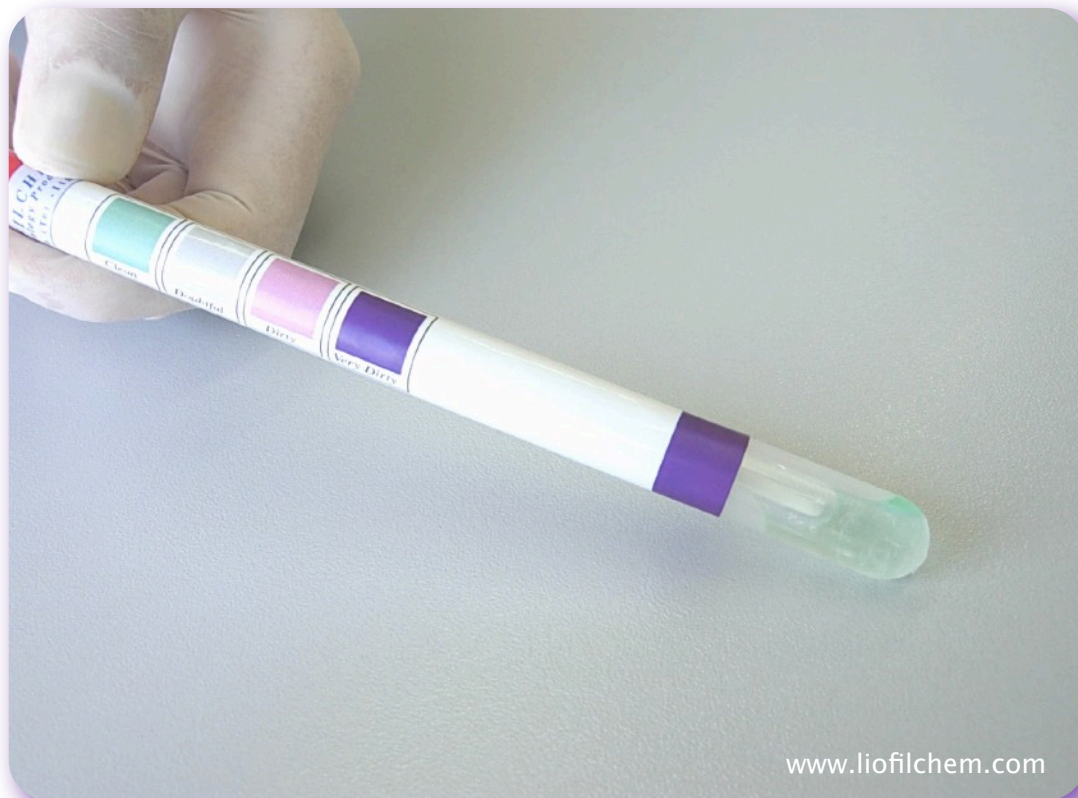

Liofilchem® and the Liofilchem company logo are registered trademarks of LIOFILCHEM s.r.l.

Swab for detecting protein, fat and sugar residues  
for the hygiene control of surfaces

Results within 10 minutes

Easy color change  
interpretation

Minimal training  
required



**Liofilchem srl**

Zona Industriale, 64026 Roseto degli Abruzzi (TE), Italy  
Tel +39 0858930745  
www.liofilchem.com  
Headquarter, Manufacturing Site, International distribution

**Liofilchem, Inc.**










465 Waverley Oaks Rd. Suite 317  
Waltham, MA 02452, USA  
Tel 781-902-0312  
US Distribution Center

ref. 6552766  
Rev. 4 - 5/12/2019

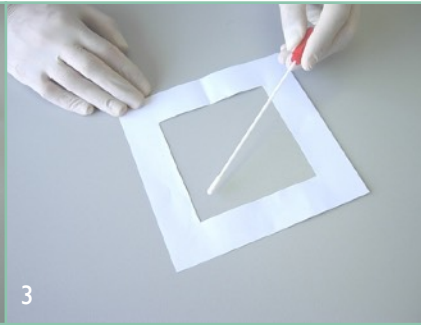

# Test procedure

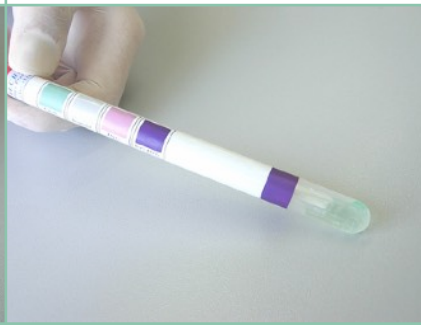Liofilchem® and the Liofilchem company logo are registered trademarks of LIOFILCHEM s.r.l.

 <p>1</p>	 <p>2</p>	 <p>3</p>
<p>The product consists of a plastic tube containing a swab, a vial (reagent A) and a disc (reagent B).</p>	<p>Take the swab contained in the tube and wet the tip with 3-4 drops of physiological solution or sterile distilled water.</p>	<p>Streak the swab on a small plot of surface measuring 10 cm x 10 cm horizontally and vertically.</p>
 <p>4</p>	 <p>5</p>	
<p>Insert the swab into the tube and push the tube in order to lower the swab tip close to the disc.</p>	<p>Break the glass vial contained in the tube with a pincer.</p>	<p>Clean surface, or dirty surface after 2 minutes reaction.</p>
		
<p>Dirty surface after 4 minutes reaction.</p>	<p>Dirty surface after 7 minutes reaction.</p>	<p>Dirty surface after 10 minutes reaction.</p>

## Clean Test

Ref. 86001

Content of the package: 50 tests



**Liofilchem srl**

Zona Industriale, 64026 Roseto degli Abruzzi (TE), Italy  
Tel +39 0858930745  
www.liofilchem.com  
Headquarter, Manufacturing Site, International distribution

**Liofilchem, Inc.**

465 Waverley Oaks Rd. Suite 317  
Waltham, MA 02452, USA  
Tel 781-902-0312  
US Distribution Center

ref. 6552766  
Rev. 4 - 5/12/2019

